

Unity Help Sheet


Renaming SpectraStar Computer for Company Networks

Some SpectraStar instruments come from the factory with a default computer name. The instructions below show how to change the name of the computer to allow for multiple SpectraStar instruments on a network or to comply with company naming policies.


1. Open **File Explorer** from the task bar.




2. Right click on **Computer** -> **Properties**.

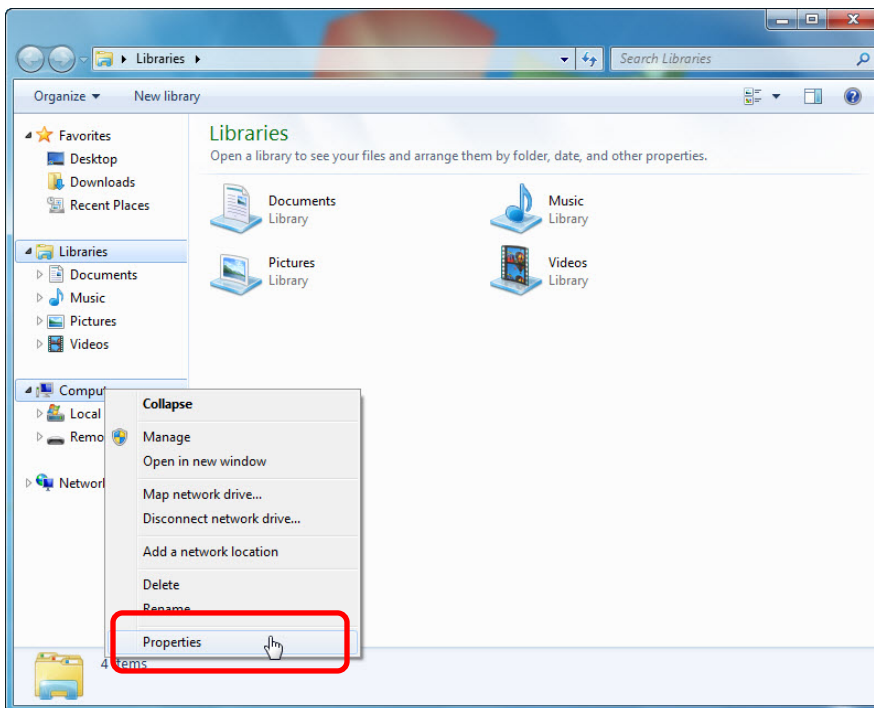


3. Press **Change Settings** on the screen below:

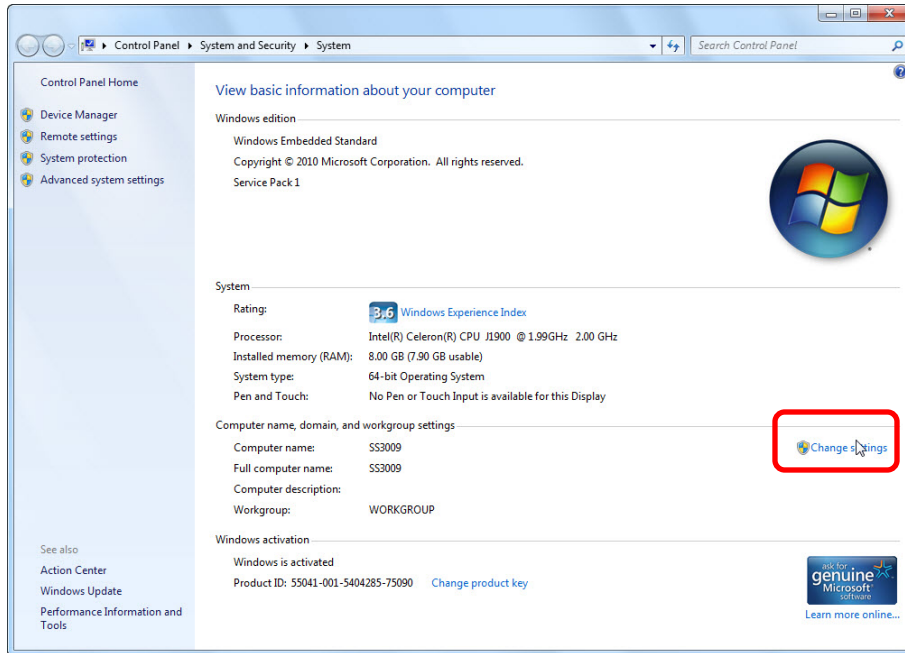


4. If desired, enter a **computer description** (not required).

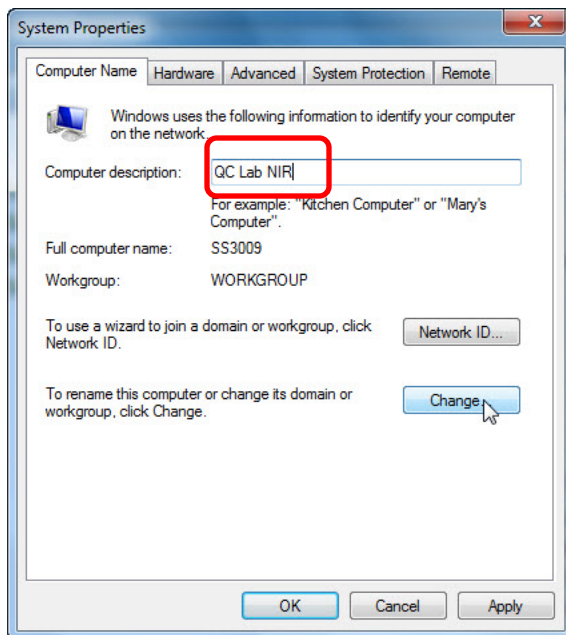



5. In the screen below, enter a **Computer name**. This name cannot contain spaces or special characters. Press **OK**.

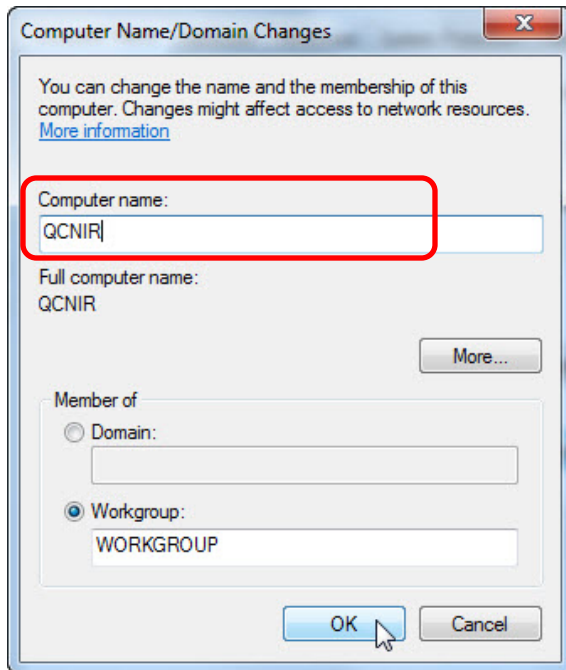

6. You will get the message below to restart the computer. Press **OK**.

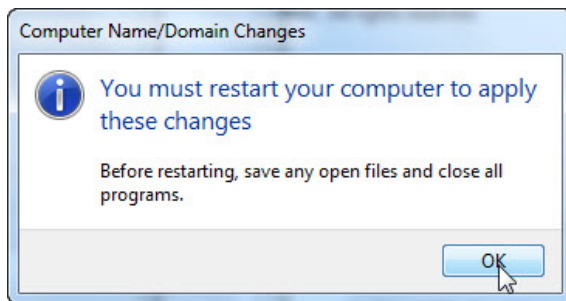

7. Press **Apply** at the next screen, followed by **Restart Now**.

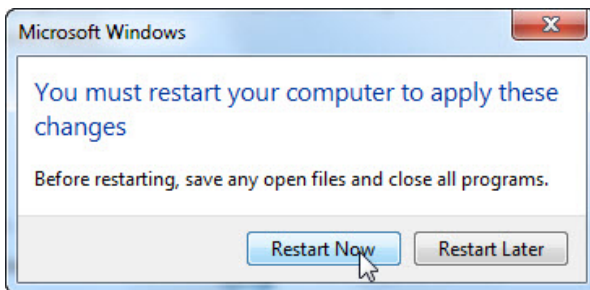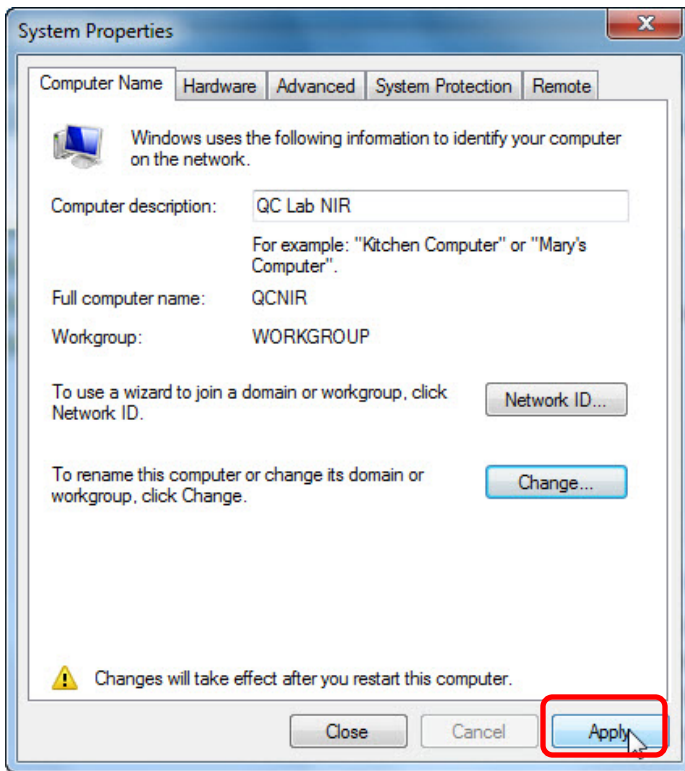

The computer name will now be changed.