


Unity Help Sheet

Updating Language Folder Permissions for Language Switching in Demo Mode


If UScan is installed in Demo mode, launching the program with the UScan Languages icon and selecting a new language can produce this error:

To solve this issue, the Windows privileges on the languages directory needs to be updated. Please follow the instructions below to update the directory rights.


1. Open **File Explorer** and navigate to **C:\Program files\Unity Scientific\UScan** directory.

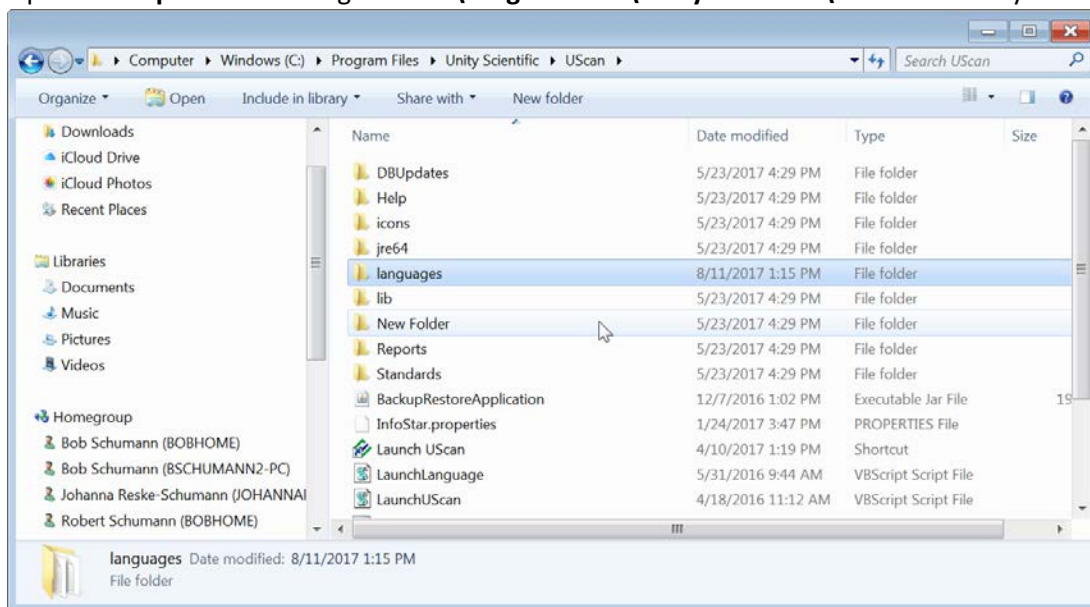


2. Right click on the languages folder -> **Properties**

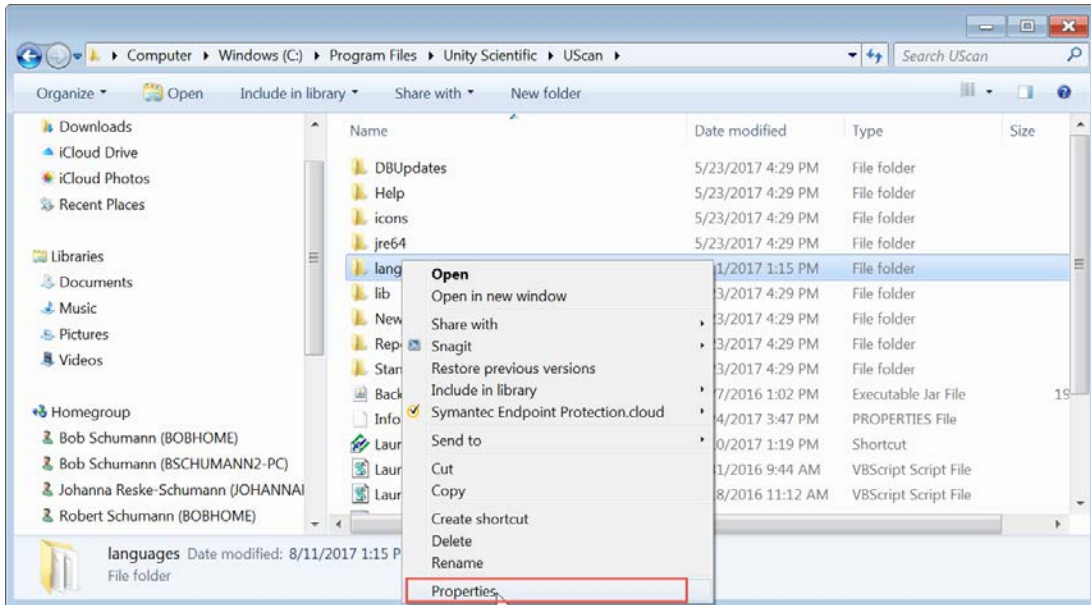

3. Click to select the **Security** tab, then **Edit** for the permissions.

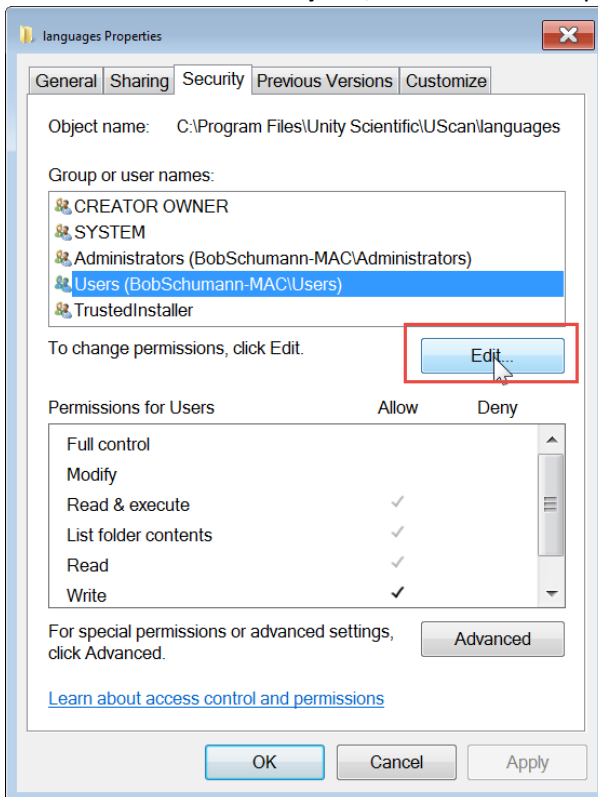

4. Click the **Users** group from the top list, then place a check mark in the **Allow** column for **Full control** and **Modify**. Press OK to close the dialog.

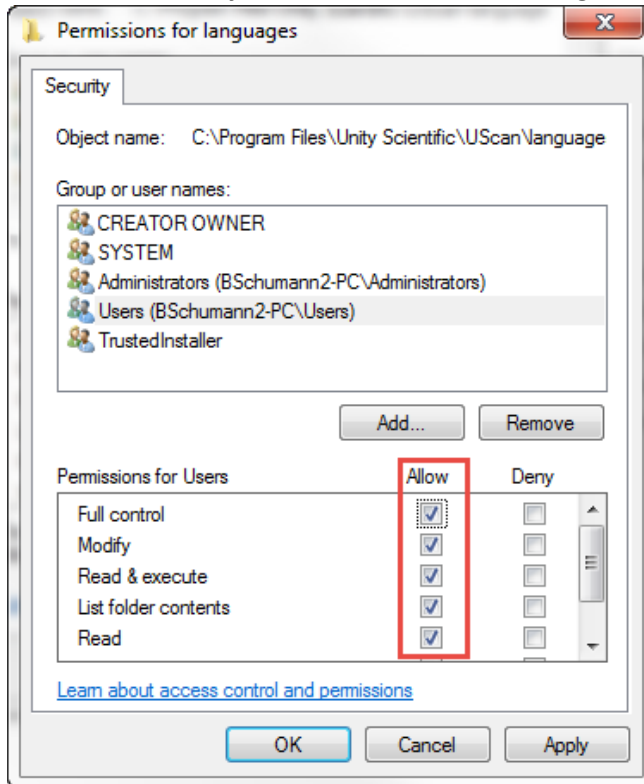

UScan should now allow you to change the language settings.